

1 CHANGING THE ZERO TRACKING TO ENABLE

Press the **[Unit]** key to view and change parameter settings when the scale is counting down from POWER UP.

The balance has 5 parameters that can be set by the user.

FUNCTION	SECTION	DESCRIPTION
UNIT F1	See section 5.1	Sets the units to be used g / N / oz
EL F2	See section 5.2	Sets the backlight AUTO EL : backlight automatically turns on when weight is applied to the top pan OFF EL : backlight always off ON EL: backlight always on
OFF F3	See section 5.3	Sets the auto power-off parameter
SETUP F4	See section 5.4	User set up parameters
tECH F5	See section 5.5	Technical parameters setting mode / factory setting

5.4 Set Up

Users can select various settings that may be suitable for certain applications. Press **[Zero/Tare]** to change a value and **[Acc]** to save changes and exit currently selected mode.

- FIL controls the speed of response of the weighing, it can be set from 0-6, default is 3. Press **[Zero/Tare]** to change, **[Acc]** is to save changes and exit this mode.
- ZEO controls the zero tracking, it can be set from 0-6, default is 3.
- ZTR controls the sensitivity, it can be set from 0-6, default is 4.
- STA controls the speed of stability indication, it can be set from 0-6, default is 2

