# MiniStirrer SM5

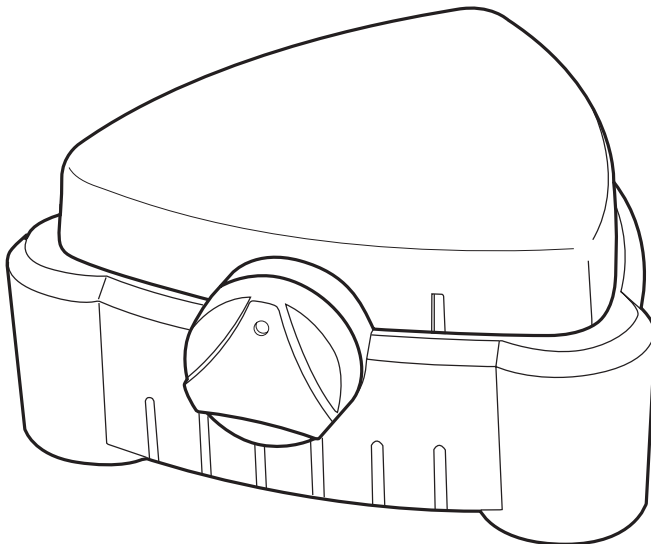
Instructions for use

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# MiniStirrer

## SM5

### Instructions for use

#### Introduction

Thank you for purchasing this piece of Stuart equipment. To get the best performance from the equipment and for your own safety, please read these instructions carefully before use. Before discarding the packaging check that all parts are present and correct.

This equipment is designed to operate under the following conditions:

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature range +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80% for temperatures up to 31°C decreasing linearly to 50% relative humidity at 40°C
- ❖ Mains supply fluctuation not exceeding 10%
- ❖ Over-voltage category II IEC60364-4-443
- ❖ Pollution degree 2

If the equipment is not used in the manner described in this manual and with accessories other than those recommended by Stuart, the protection provided may be impaired.

#### Electrical Installation

**THIS EQUIPMENT MUST BE EARTHED**

**Before connection please ensure that the line supply corresponds to that shown on the rating plate.**

**Model Supply requirements**  
SM5 100-230V-AC, 50/60 Hz

All models are supplied with a mains supply lead fitted with a 12 Volt plug for connection to the unit.

**NOTE:** Refer to the equipment's rating plate to ensure that the plug and fusing are suitable for the voltage and wattage stated.

**IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN**

#### Safety Advice Before Use

- ❖ The unit is not spark, flame or explosion proof and has not been designed for use in hazardous areas in terms of BSEN 60079-14:1997. Keep flammable, low flash point substances away from the apparatus.
- ❖ This equipment will generate magnetic fields. Keep all magnetic data devices (e.g. credit cards) away from the stirrer unit. The effects of the magnetic field have to be taken into account with respect to cardiac pacemakers etc.
- ❖ **DO NOT** immerse in water or fluids.
- ❖ Avoid surface contact with aggressive media.

#### General Description

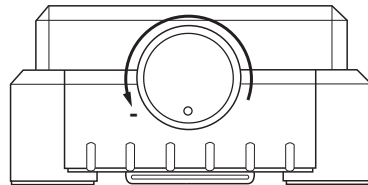
The MiniStirrer is designed for stirring fluids in a flat bottomed glass vessel using a magnetic stir bar. It is available in three fascia designs: SM5/STUART, SM5/SWIRL and SM5/BIBBY. It can stir up to 1 litre of fluid in the range 350 to 2000rpm.

#### Installation


Locate the MiniStirrer on a solid work surface with 300mm clearance from other equipment. Ensure the power disconnect device is accessible.

#### Operation


1. Rotate the control knob **anti-clockwise** to the fully off position.

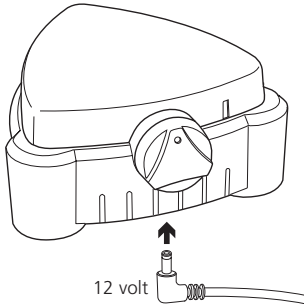
2. Insert the power supply 12 Volt plug into the base socket of the MiniStirrer.

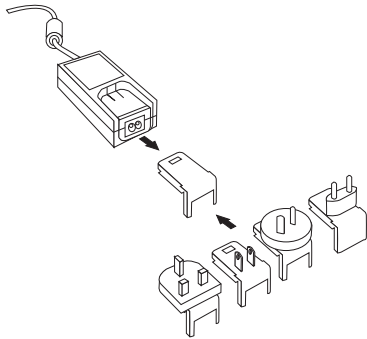
3. Select the correct plug attachment and attach to the power supply unit as shown below. Plug the power supply plug into the mains supply socket.

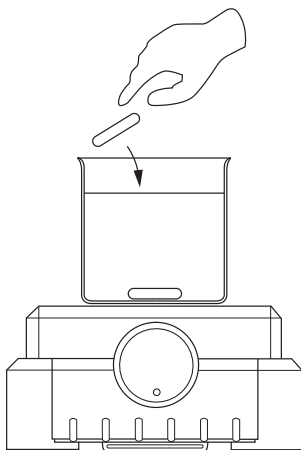4. Place a loaded beaker containing a magnetic stir bar on the MiniStirrer platform.

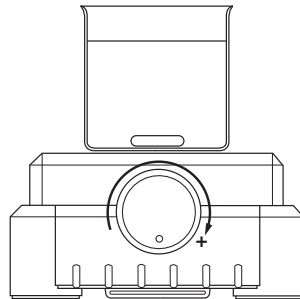5. Rotate the control knob **clockwise** to adjust the stirrer speed between 350-2000rpm.

## Maintenance & Servicing

The MiniStirrer **is not** serviceable. Routine maintenance should include inspection of the power supply unit cord and stirrer base for damage. Periodically clean the instrument using a dry cloth.

For any technical enquiries regarding this product, please contact the Technical Support Department:  
E-mail: [stuarthelp@bibby-scientific.com](mailto:stuarthelp@bibby-scientific.com)  
Tel: +44 (0)1785 810433

## Warranty

Stuart warrants this equipment to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of **three (3)** years. In the event of a justified claim, Stuart will replace any defective component or replace the unit free of charge. This warranty does NOT apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage caused by installation, adaptation, modification, fitting of non-approved parts or repair by unauthorised personnel.

Bibby Scientific Ltd.  
Beacon Road, Stone, Staffordshire  
ST15 0SA  
United Kingdom  
Tel: +44 (0) 1785 812121  
Fax: +44 (0) 1785 813748  
E-mail: [equipment@bibby-scientific.com](mailto:equipment@bibby-scientific.com)  
[www.stuart-equipment.com](http://www.stuart-equipment.com)

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# Notes

# Notes

These products meet the relevant EC harmonised standards for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in their vicinity will meet these standards

and we cannot guarantee that interference will not occur in practice. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advice before use, please contact the Technical Service Department of Bibby Scientific Ltd.

## Declaration of Conformity

### Laboratory Mini Stirrer, Model SM5

These products comply with the requirements of the EU Directives listed below:

**2004/108/EC**      **EMC Directive.**  
**2006/95/EC**      **Low voltage Directive (LVD)**

Compliance with the requirements of these Directives is claimed by meeting the following standards:

**EN 61326-1:2006 (Electrical Equipment for Measurement, Control and Laboratory use).**  
**EN 61010-2-051 : 2003 (Safety requirements for electrical equipment for measurement, control and laboratory use. Particular requirements for laboratory equipment for mixing and stirring)**  
**FCC 47CFR Part 15B Class B**

CE mark affixed '12

Signed:  (Mr S. Marriott)

Date: 1/2/12

Authority: Technical Director  
Bibby Scientific Ltd

Bibby Scientific Ltd - Stone - Staffs - ST15 0SA - UK  
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

**UK**

Bibby Scientific Ltd.  
Beacon Road, Stone  
Staffordshire ST15 0SA  
United Kingdom  
Tel: +44 (0)1785 812121  
Fax: +44 (0)1785 810405  
e-mail: [sales@bibby-scientific.com](mailto:sales@bibby-scientific.com)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)

**France**

Bibby Scientific France SAS  
ZI du Rocher Vert - BP 79  
77793 Nemours Cedex  
France  
Tel: +33 1 64 45 13 13  
Fax: +33 1 64 45 13 00  
e-mail: [bsf@bibby-scientific.fr](mailto:bsf@bibby-scientific.fr)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)

**North & South America**

Bibby Scientific US Inc.  
t/a Techne Inc.  
3 Terri Lane, Suite 10  
Burlington, NJ 08016 USA  
Toll Free (in NA): 800-225-9243  
Tel: +1 609 589 2560  
Fax: +1 609-589-2571  
e-mail: [labproducts@techneusa.com](mailto:labproducts@techneusa.com)  
[www.techneusa.com](http://www.techneusa.com)

**Italy**

Bibby Scientific Italia Srl  
Via Alcide de Gasperi 56  
20077 Riozzo di Cerro al Lambro  
Milano Italia  
Tel: +39 (0)2 98230679  
Fax: +39 (0)2 98230211  
e-mail: [marketing@bibby-scientific.it](mailto:marketing@bibby-scientific.it)  
[www.bibby-scientific.it](http://www.bibby-scientific.it)

**Middle East**

Bibby Scientific Middle East Ltd.  
PO Box 27842, Engomi 2433  
Nicosia  
Cyprus  
Tel: +357 22 660 423  
Fax: +357 22 660 424  
e-mail: [sales@bibby-scientificme.com](mailto:sales@bibby-scientificme.com)

**Asia**

Bibby Scientific - Singapore  
Prudential Tower, Level 26  
30 Cecil Street  
Singapore 049712  
Tel: +65 6631 2976  
Fax: +44 (0)1785 810405  
e-mail: [info@bibby-scientific.com](mailto:info@bibby-scientific.com)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)